

### **Transferring Liquids Using a Calibrated Pipette**

Place the bulb on the end of the pipette. Place the tip of the pipette in the liquid and squeeze slightly so the liquid will come up into the glass pipette slightly above the level you need, but not into the bulb. Remove the bulb and replace with your index finger. Release the liquid you do not need by removing your index finger slightly until the level of the liquid is at the volume mark that you wish to transfer. Hold the pipette over the vessel to which it is to be transferred. Release your index finger.

Wash the pipette after use by squeezing water into pipette (not the bulb) and releasing it into the sink. Repeat several times.